GCWPS Cartridge Filters

High Capacity PES Membrane





GCWPS Membrane Filter Cartridges are designed for general service wherever a membrane filter is required. Designed to hold the maximum amount of filter media that can be effectively used in a cartridge, GCWPS filters lower the cost of filtration. GCWPS cartridges are flushed with high purity water to remove manufacturing debris.

Construction Materials

Filtration Media	High Capacity Polyethersulfone (PES) Membrane		
Media Support	Polypropylene		
End Caps	Polypropylene		
Center Core	Polypropylene		
Outer Support Cage	Polypropylene		
Sealing Method	Thermal Bonding		
O-rings	Buna, Viton® (or FKM), EP, Silicone, FEP Encapsulated Silicone, FEP Encapsulated Viton (or FKM)		

Dimensions

Length	5 to 40 in. (12.7 to 101.6 cm) nominal
Outside Diameter	2.75 in. (7.0 cm) nominal
Filtration Area	6.0 ft² (0.56 m²) per 10 in. length

Applications

- Process Water
- **♦** Chemicals
- DI Water
- Cosmetics
- Inks and Dyes

Integrity Test Information

Representative samples from each manufacturing lot are tested for integrity to ensure consistent performance.

Maximum Operating Parameters

Differential Pressure • Forward	50 psid (3.4 bard) at 20 °C (68 °F)		
• Reverse	40 psid (2.7 bard) at 20 °C (68 °F)		
Operating Temperature	82 °C (180 °F) at 10 psid (0.69 bard) in water		
Recommended Changeout Pressure	35 psid (2.4 bard)		

Sanitization/Sterilization

Filtered Hot Water	90 °C (194 °F), 30 minutes, multiple cycles, max 3 psid forward flow
Autoclave	121 °C (250 °F), 30 min, multiple cycles
In-line Steam	135 °C (275 °F), 30 min, multiple cycles

For all elevated temperature procedures above, a stainless steel support ring is required.

Chemical Sanitization

Performed using industry standard concentrations of hydrogen peroxide, paracetic acid, sodium hypochlorite and other selected chemicals.

Total Performance

Critical Process Filtration, Inc. is a vertically integrated manufacturer of filtration products to industries in which filtration is considered a critical part of the manufacturing process. We supply a complete line of products and services to help you cost effectively satisfy all your filtration requirements from a single source.

FDA and EC Compliance

All Critical Process Filtration filters are designed to meet the FDA requirements for processing food and beverage products. The materials used to construct GCWPS filters are listed by the FDA as appropriate for use in articles intended for repeated food contact as specified in Title 21 CFR sections 174.5, 177.1500, 177.1520, 177.1630, 177.2440 and 177.2600 as appropriate. Membrane filters comply with Title 21 CFR sections 210.3 (b)(6) and 211.72, for nonfiber releasing filters. All materials used to make the filters are listed in European Commission Regulation EU/10/2011, Annex 1.

Quality Assurance and Standards

Our goal is to ensure our customers the greatest possible value for their filtration dollar. Our state of the art manufacturing facility and quality management system both meet ISO 9001:2008 standards. Each operation from assembly and test to cleaning, drying, and packaging is done in appropriately rated clean rooms. A sophisticated MRP system collects and processes real time data from manufacturing centers and inspection points. This allows variable and attribute data to be quickly and easily analyzed driving constant improvements in both quality and cost.

Extractables

GCWPS filters generally exhibit low levels of non-volatile residues.

Flow Rate


The Typical Flow Rates table represents typical water flow at a 1 psid (69 mbard) pressure differential across a single 10 in. cartridge element. The test fluid is water at ambient temperature. Extrapolation for housings with multiple elements and higher pressure drops is acceptable, but as flows increase the pressure drop of the housing becomes more apparent.

Typical Flow Rates

Pore Size	0.22 μm	0.45 μm	0.65 μm	1.2 μm
GPM	1.1	1.4	4.5	7.0
LPM	4.16	5.30	17.03	26.50

Ordering Information

Cartridge order numbers have several variables from pore size to end cap type. For example, General Service Grade High Capacity PES Membrane, 0.22 Micron Rating, No SS Support Ring, 20"Length, Silicone O-Rings, 2-222/Flat End Cap Configuration = GCWPS-20N00002S5.

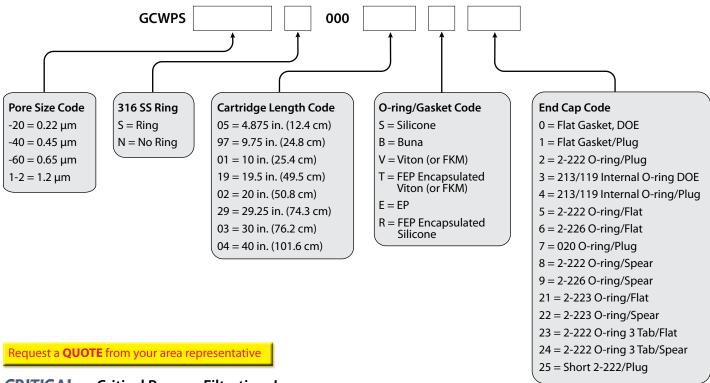

CRITICAL PROCESS I I I I I FILTRATION, INC.

Critical Process Filtration, Inc.

One Chestnut Street • Nashua, NH 03060 Tel: 603.880.4420 • Fax: 603.880.4536

criticalprocess.com • sales@criticalprocess.com

The information contained herein is subject to change without notice.

The Critical Process Filtration logo is a trademark of Critical Process Filtration, Inc.

Viton is a trademark of DuPont Performance Elastomers L.L.C.

© 2012-2015 Critical Process Filtration, Inc. • All Rights Reserved • Data Sheet GCWPSDS1211 Rev-